



Control D Level 2 Bio

for Cedex Bio and Bio HT Analyzers

REF 07 368 186 001

6 × 1.0 mL

Control Code: COND2, 308, 07-1144-6

Version 14
Content version: December 2019

Store at +2 to +8°C

Intended Use

This product is intended for quality control by monitoring accuracy and precision of quantitative assays using Cedex Bio and Bio HT Analyzers.

Use this control as described in the Instructions for Use for the respective assay.

Storage and Stability

Store at +2 to +8°C.

Stable at +2 to +8°C until the expiry date printed on the label when stored unopened in original vials and kept free of contamination.

Open vial is stable for 1 day at +15 to +25°C.

Procedure

The control is ready for use. Mix carefully before use.

Avoid formation of foam. Air bubbles can be removed by centrifugation in a microcentrifuge for 10 to 20 seconds.

Changes to Previous Version

- New lot-specific information.

Trademarks

CEDEX is a trademark of Roche.

Online Technical Support

Please visit our Online Technical Support site for additional information about this product: www.technical-support.roche.com

Lot-Specific Data

Lot No.: 46341900

Expiry Date: Dec 2020

Assigned concentration values:

Analyte	Value	1 s	Unit
Acetate	18.738	0.937	mmol/L
	1106.3	55.3	mg/L
Ala-Gln	4.000	0.200	mmol/L
	868.9	43.4	mg/L
Formate	8.000	0.400	mmol/L
	360.2	18.0	mg/L
Galactose	6.994	0.350	mmol/L
	1260.0	63.0	mg/L
Pyruvate	3.000	0.150	mmol/L
	261.1	13.1	mg/L
Sucrose	16.400	0.820	mmol/L
	5613.7	280.7	mg/L

**For more information please visit
custombiotech.roche.com**

© 2019 Roche Diagnostics GmbH. All rights reserved.

For more information, contact your local Roche representative.

Europe, Middle East, Africa,
and Latin America:
Phone +49 621 759 8580
Fax +49 621 759 6385

Asia Pacific:
Phone +65 6371 6638
Fax +65 6371 6601

Japan:
Phone +81 3 6634 1046
Fax +81 3 5479 0585

Canada:
Phone +1 450 686 7050
Fax +1 450 686 7012

United States:
Phone +1 800 428 5433,
ext. 14649 (toll-free)
Fax +1 317 521 4065

Roche Diagnostics GmbH
Sandhofer Str. 116
68305 Mannheim
Germany

